

Vacuum Manifold Instructions

1. Remove caps from enough ports to hold all of your samples
2. Place columns onto ports and make sure valves are open
3. Place samples into columns
4. Turn vacuum on until all samples have flowed through column
5. Turn vacuum off
6. Repeat steps 3-5 as needed to complete protocol
7. Once all samples are completed, replace caps over ports
8. Remove hose and take manifold to sink
9. Remove black plug and pour flow through liquid down sink
10. Replace plug securely, wipe down manifold
11. Return to counter and replace hose
12. Wipe down bench area with 70% ethanol and take all supplies back to your bench

